# NKT Photonics

• •

# Koheras HARMONIK HPH76 High-power UV & VIS fiber laser system



### A HAMAMATSU COMPANY

# Single frequency, low noise

## Ideal for cutting-edge quantum physics projects

HARMONIK HP is a range of high-power frequency-converted fiber laser systems with low linewidth and phase noise. Upon request, further linewidth and RIN reduction options are available.

### High output power and excellent beam quality

The output power of the HARMONIK HP systems depends on the wavelength, with the highest obtainable output of 10 W. The excellent beam quality,  $M^2 < 1.1$ , is suitable for cutting-edge quantum physics projects such as quantum sensing and laser cooling and trapping.

### Ideal for quantum optics

With its ultra-stable and narrow linewidth, the HARMONIK HP is ideal for applications such as quantum optics, computing and other phenomena like optical trapping, optical lattice, Bose-Einstein condensate, atom interferometer, and squeezing.

### Rack system and table top

The BOOSTIK HP system is delivered in a turn-key 19" rack system and the HARMONIK HP frequency-conversion modules are supplied as tabletop units that also fit a 19" rack-mount module.



Koheras HARMONIK HP H76

## Koheras HARMONIK HP H76

## Applications

Quantum computing **Atomic trapping and cooling** 

BENEFITS

## Benefits

### Ultra-low noise

The HARMONIK HP systems offer an ultra-low phase noise as well as low RIN. The RIN is preserved from the pump laser while the linewidth is doubled.

### Fast wavelength modulation

The systems are equipped with easy and user-friendly fast wavelength modulation. This feature is typically used to lock the laser to an external stable reference — such as an atomic transition line or interferometer — to obtain even higher wavelength stability than provided by the free-running laser.

### Thermal tuning

All Koheras fiber lasers are equipped with thermoelectric temperature controllers (TECs).


The TECs stabilize the operation of the laser and makes it insensitive to environmental temperature fluctuations. The TECs also make it possible to tune the center wavelength by changing the operating temperature of the laser.

At standard room temperature (around 20-30°C or 68-86°F) the laser can be thermally tuned to an industry-leading 500 pm.

### Free-space or fiber output

As default, the HARMONIK HP systems come with free-space output. If fiber coupling is needed, we can propose an efficient coupling via our unique polarization-maintaining single mode photonic crystal fiber that ensures high-power delivery.

Our pre-angled fiber couplers are easy and fast to align. They give a very high coupling efficiency, which does not decrease significantly over time, so it is possible to obtain a long term stable coupling.

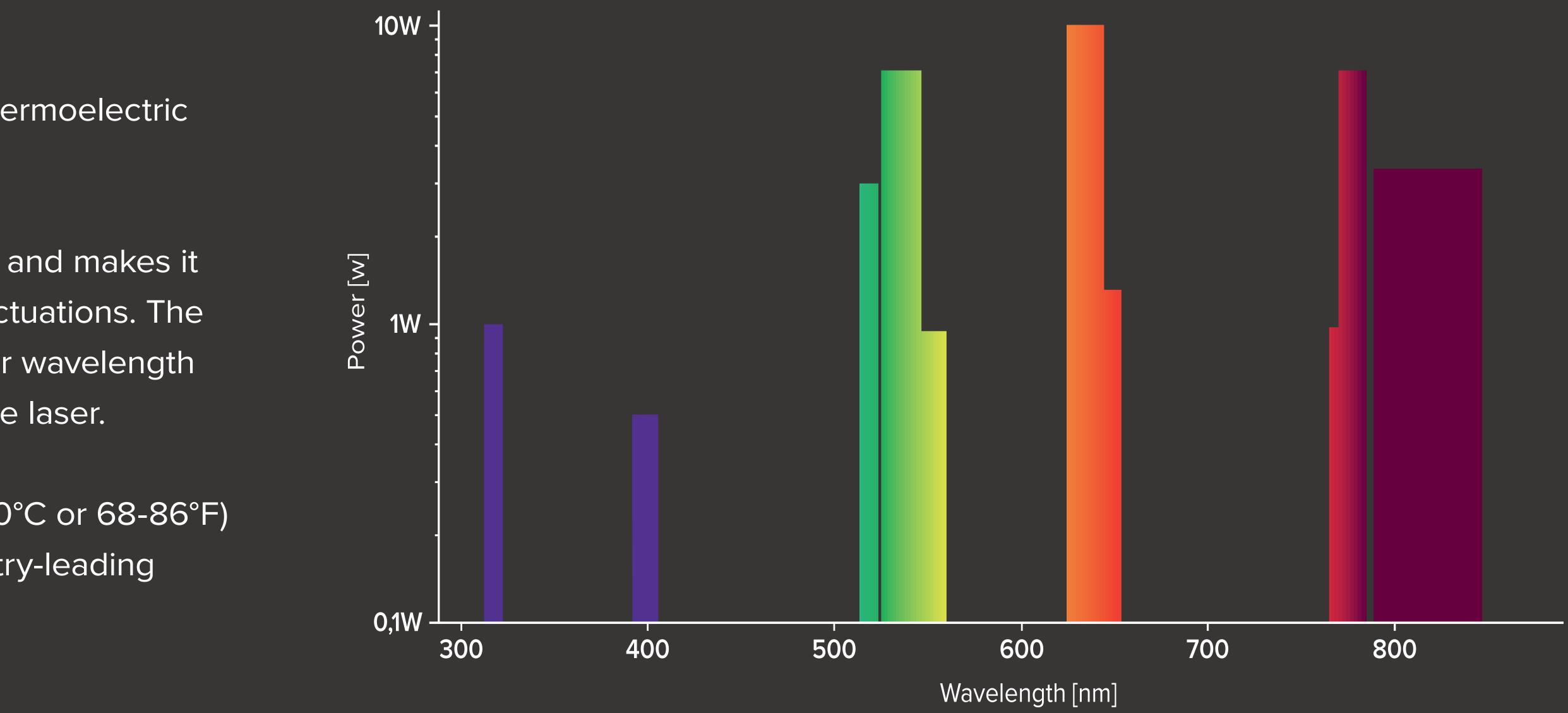

## Koheras HARMONIK HP H76

The graph shows the VIS wavelength capabilities of the HARMONIK HP.

All lasers are pumped by our low-noise fiber lasers in the NIR (not shown), allowing the lasers to be locked to frequency references at either their fundamental or converted wavelengths.

## Specifications

## Optical

Model

Center wavelength <sup>1</sup> [nm]

Standard wavelength [nm]

Laser emission

**Beam quality** 

Output power <sup>2</sup> [W]

Linewidth [kHz]

RIN peak [MHz]

RIN level [dBc/Hz]

@ peak

@ 10 MHz

Long term stability (peak-peak, 1h @25°C) <sup>3</sup> [%]

Optical S/N (50 pm res.) [dB]

Polarization / PER [dB]

Min. thermal wavelength tuning range [pm]

Total thermal wavelength tuning range [pm]

H76	Fast wavelength modulation range [GHz]	> 16
765-769	Fast wavelength modulation [kHz]	Up to 20
767	Pump unit(s) size for 19" rack system [U]	4
CW - Inherently single frequency	Number of frequency conversion modules <sup>4</sup>	1
M <sup>2</sup> < 1.1		
> 4	1) Center wavelength is selectable within the specified range. 2) Shot noise-limited > 5 Mhz. 3) After a 10-20 minute warm-up.	
< 0.4	4) Delivered as tabletop module(s).	
≈ 0.7		
< -97		
< -132		
< 10		
> 70		
> 20		
± 175		
500		

Koheras HARMONIK HP H76

## Koheras HARMONIK HP H76

## Service and warranty extensions

The Koheras warranty and service package ensures trouble free operation of your Koheras laser. The Standard Package gives you a two year warranty extension plus remote diagnostics of key laser parameters through a remote connection to the system.

Our Premium Package adds a guarantee that we always stock a laser with your specifications - ready to ship - should you need it.

## Specifications

## Electrical/Mechanical/Environmental

### HARMONIK HP H76<sup>1</sup>

Power supply requirements [VAC, Hz]

**Digital interface** 

Monitor output

**Dimensions (WxHxL) [mm]** 

Weight [kg]

1) For the BOOSTIK HP system, please refer to the BOOSTIK HP datasheet.

100-240 VAC, 50-60 Hz

Ethernet 10/100

Standard: Free-space

Optional: 2.5m high power, single mode fiber delivery FC/APC

173 x 62.5 x 240

6

Koheras HARMONIK HP H76


## Koheras HARMONIK HP H76


## Reliability


The Koheras range of single frequency fiber lasers is based on telecom-grade fiber components and built to last thousands of hours with no service or maintenance.


With several thousand lasers installed in environments varying from fully climate controlled national standards laboratories to the demanding environment on oil rigs and submarines, the Koheras line is the most robust singlefrequency laser range on the market with an unmatched reliability track record.

## Technical Drawings

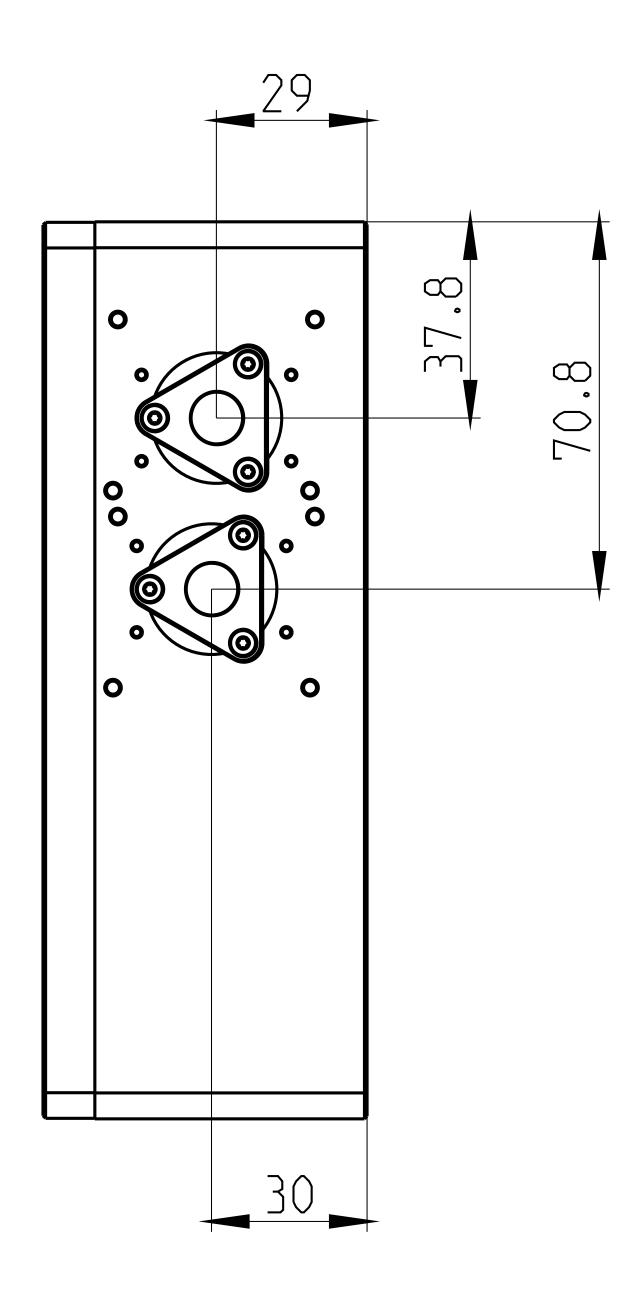

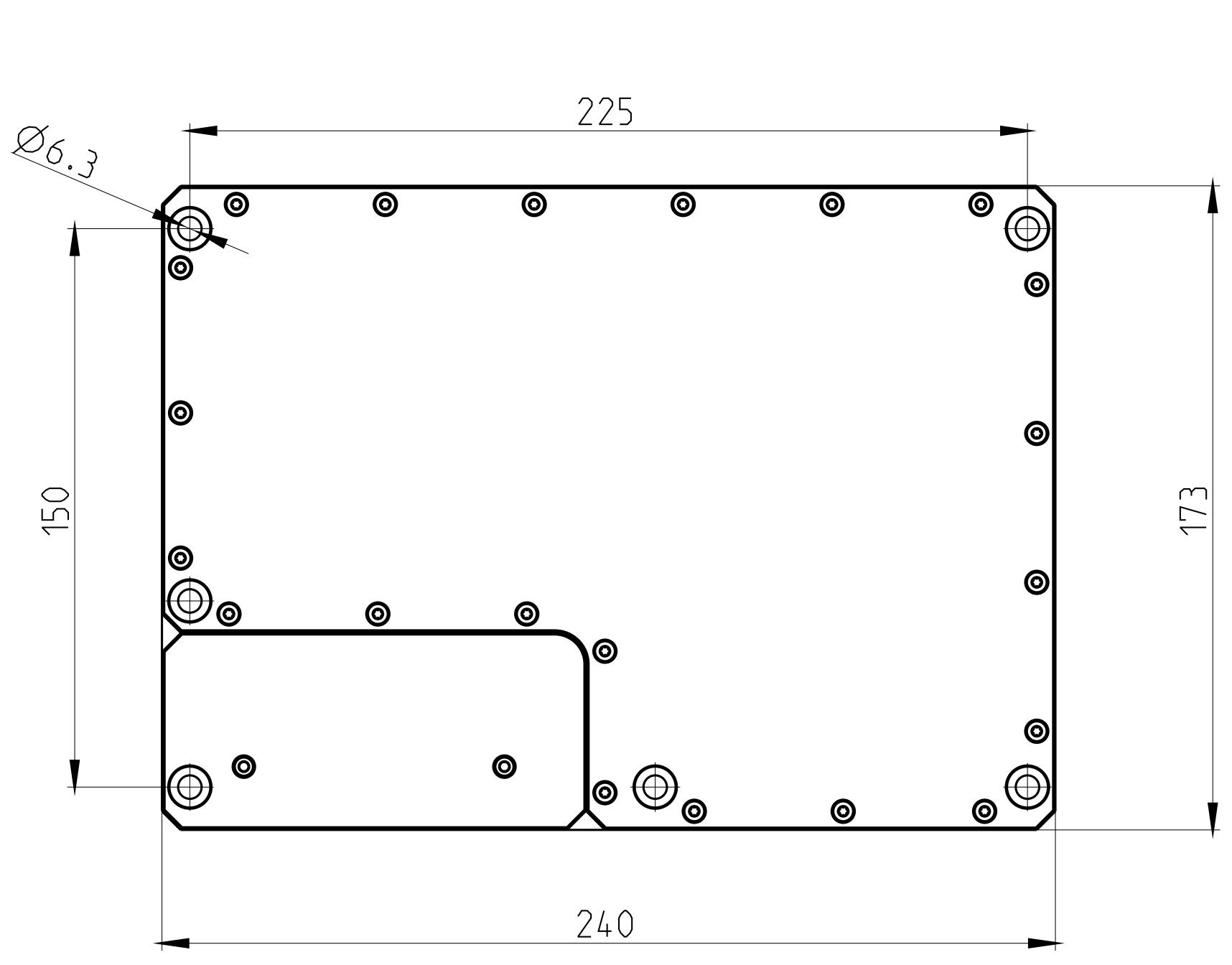

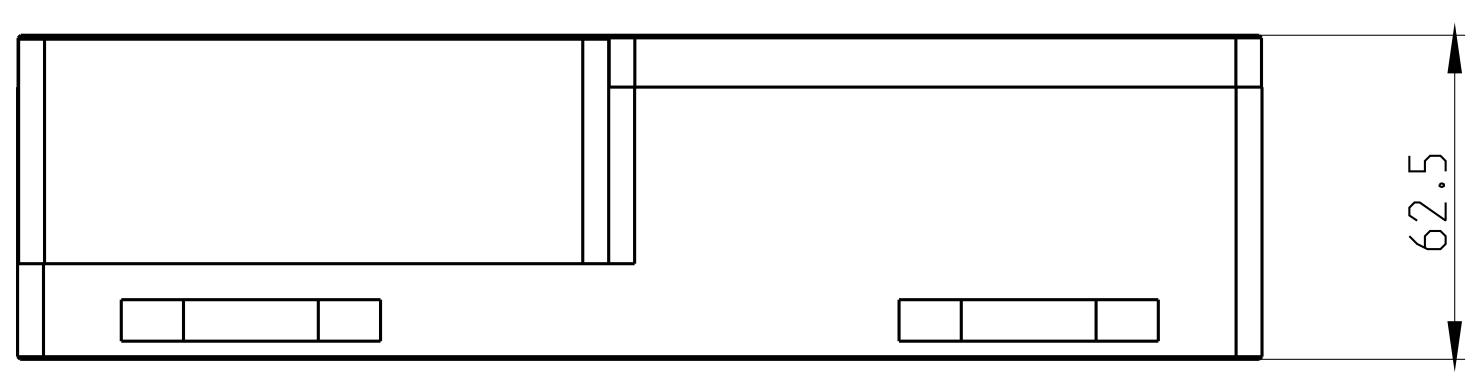

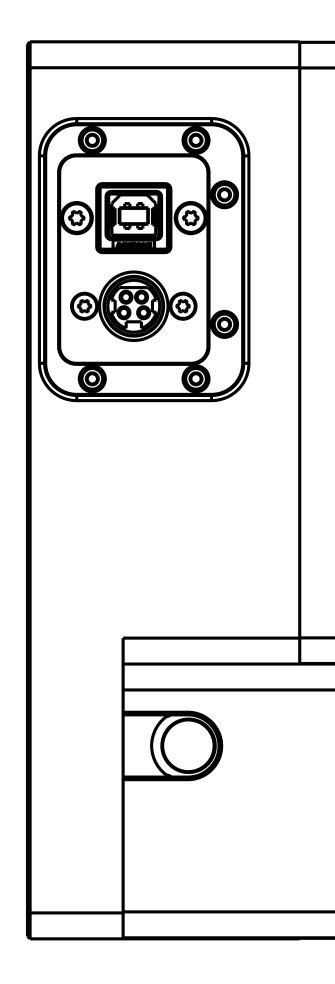

## Koheras HARMONIK HP H76

All NKT Photonics products are produced under our quality management system certified in accordance with the ISO 9001:2015 standard.









### A HAMAMATSU COMPANY